

# CELLULAR STAINING PROTOCOL


## N-Propanoyl-BODIPY-monosialoganglioside GM<sub>1</sub> (NH<sub>4</sub><sup>+</sup> salt)

Catalog number 2060

### Staining SH-SY5Y Cells with C3-BODIPY-GM<sub>1</sub>

1. Dissolve C3:0-BODIPY-GM<sub>1</sub> (100 µg) in 64 µl DMSO:methanol:water (2:1:0.2) and aliquot. Store aliquots at -20°C until use.
2. Plate cells at density of 10,000/well in a 96-well plate in media with 10% FBS and allow to attach overnight.
3. The following day, remove media and replace with 2.5-10 µM C3:0-BODIPY-GM<sub>1</sub> in phenol red-free media. Return cells to incubator for 30 minutes.
4. Remove media and replace with fresh phenol red-free media. Image using a fluorescence microscope with a FITC/GFP laser.

Note that complexing fluorescent lipids with bovine serum albumin (BSA) facilitates cell labeling by eliminating the need for organic solvents to dissolve the lipophilic probe. A BSA-complexed probe can be directly dissolved in water. However, gangliosides are unique lipids that are soluble in aqueous systems.

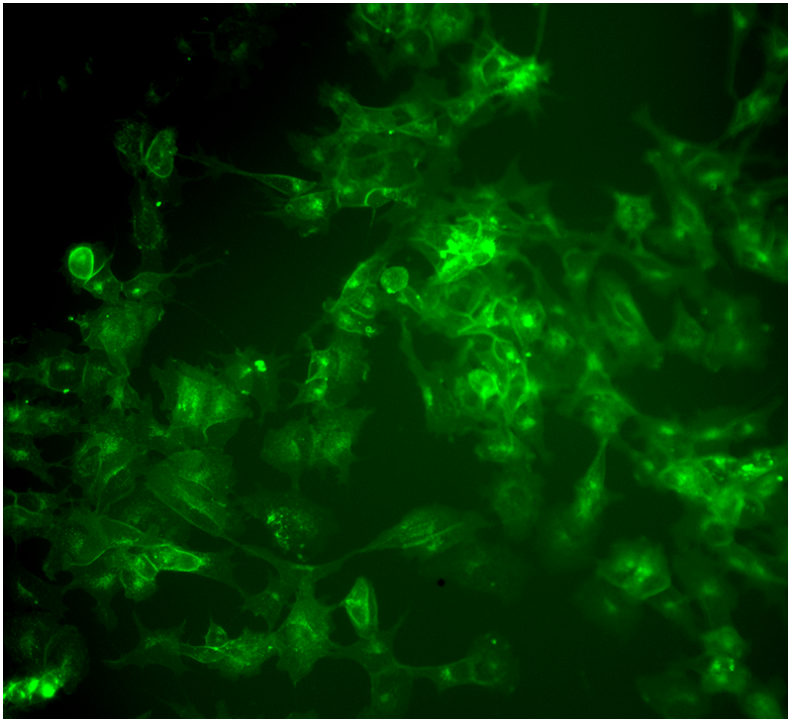

SH-SY5Y neuroblastoma cells stained with 10 µM C3-BODIPY GM<sub>1</sub>.

This product is to be used for research only. It is not intended for drug or diagnostic use, human consumption, or to be used in food or food additives. Matreya assumes no liability for any use of this product by the end user. We believe the information, offered in good faith, is accurate.